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Stock List for Urgent service

Tag Name	Name	AL Part No	Type	Supplier	Description	Application model	Reco. stock	Priority
FS615/635 for Seal	Flowswitch		SMR12GGXFRKG/U S-100 25+ ADOAH040VAS0005 E04 M12 +E40199	ifm efector	Type: Flow meter SM6000 ,End Connection:ISO BSP1/2" Outs. Thr., Material Product Wetted Part:EN 1.4404(SS316L), Measuring Range :(0.1-25)L/Min, Cable Entry:M12 Connector, Power supply: 30DC ,Output:0-10V +4-20 mA ,Angled connector EVT004 with 5 mt.cable ,Adapter REDADAPT G1/2"- R1/2"	Brew250/Brew2000/Brew 3000 /CR250/PX810(Bactofuge) /Clara H series	1	1
PT 375 for OP water	Pressure Transmitter	569089-01	MBS 3000-111-AB04	Danfoss	Type: Pressure Transmitter MBS 3000, End Connection: ISO 228 BSP 1/4" Welded,Material Product Wetted part: Stainless Steel, Measuring Range:0/2.5,Cable Entry: Power Supply 28VDC	Clara200/CR250 /Clara series	1	1
ST	RPM sensor for HSS	552042-01	Inductive Sensor Namur	PEPPERL+FUCHS	PEPPERL+FUCHS NCN4-12GM35-NO 10 M	HSS	1	1
ST	RPM sensor for HSS	552042-02	Inductive Sensor Namur	PEPPERL+FUCHS	PEPPERL+FUCHS NCN4-12GM35-NO 2 M	Clara20	1	
ST	RPM sensor for Old Decanter	61208903-01	Inductive Sensor Namur	PEPPERL+FUCHS	PEPPERL+FUCHS 2 wire	Old Decanter	1	
ST	RPM Sensor for 2 touch	61211273-30	Inductive Sensor Namur	PEPPERL+FUCHS	PEPPERL+FUCHS 2wire	Old 2 -touch	1	1
ST	RPM Sensor for 2 touch plus	61244548-01	Inductive Sensor Namur	ifm efector	IGT 219 3 wire	Foodec 300 /P2-705	1	
Inlet seal	CF100	585712-81	Inlet seal	AlfaLaval	KOL/SIC-KOL/SIC FDA	CF100 (6 units)	1	1
YT	Vibration Sensor	61244012-01	VKV022	ifm efector	VKV022 Vibration Velocity RMS	Old Decanter & Old 2 touch	1	1

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Click!
Pictures to
view files

Q&A for Biotech/Pharma HSS

Subject		Document No.	Rev. no.
Operator training of Biotech/Pharma HSS			
Issued by	Department	Date	Approved by
DaeKi Hong	Service	2022/10/11	
Classification			

Q&A Biotech /Pharma HSS

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Factor	Before	After	Remark
Flow rate			
Discharge Interval Or Discharge Type			
Back Pressure			
Flow control mode			
Solid Volume			
Temp			
IP Conveter Function			
Water pressure			
Discharge Size			

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Factor	Before	After	Remark
Vibration Level			
Discharge Time & Size			
Feed condition			
Motor Amp			
Noise Position			
Exchange IS Kit			
Exchange MS Kit			
Exchange Oil			
Safety water			

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